

## Attune NxT 2 Configuration

	Mirror	Filter	Fluorochrome Examples
<b>488</b>	503LP	530/30	FITC, GFP, AF488
50mw	555LP	695/40	PerCP-Cy5.5
<b>637</b>	654LP	670/14	APC
100mw	690LP	720/30	AF700
	740LP	780/60	APC-Cy7
<b>405</b>	417LP	450/40	Brilliant Violet 421, eFluor 450, Pacific Blue, DAPI
50 mw	495LP	525/50	eFluor 506, Brilliant Violet 510, Pacific Green
	555LP	610/20	Brilliant Violet 605, Pacific Orange
	635LP	660/20	Super Bright 645, Brilliant Violet 650
	680LP	710/50	Super Bright 702, Brilliant Violet 711
	740LP	780/60	Brilliant Violet 786
<b>561</b>	577LP	585/16	PE, dTomato, DsRed
50 mw	600LP	620/15	PE-Texas Red, AF594, mCherry
	650LP	780/60	PE-Cy7

Detector	Starting Voltages (Using Beads and Lymphocytes)
BL1	350
BL2	365
RL1	420
RL2	440
RL3	425
VL1	300
VL2	330
VL3	375
VL4	365
VL5	390
VL6	400
YL1	380
YL2	420
YL3	410

Update: 09-25-2025