



ATTUNE USER GUIDE

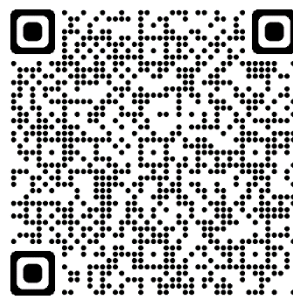

ILABS KIOSK

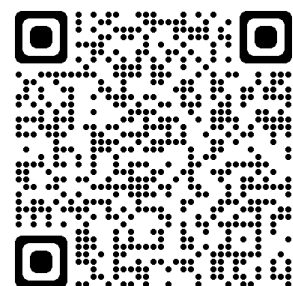

If issues aren't resolved,
report to Staff

Instrument Issue Reporting



QUICK START

Start Up (First User)

Start Kiosk session

Power cycle instrument. Make sure to turn Autosampler on FIRST then Instrument

Log on¹ and Run Start up in Instrument Tab

Run PT test

Experiment Set up

Create New Experiment
(Template or New)

Verify *Instrument Settings*

Check controls and adjust settings accordingly

Compensation

Use "Unstained" Option if you do not have negative populations in tubes

Use "Negative Gate" if negative populations exist in ALL samples

Record compensation controls

Acquisition

Create Groups/Wells/Tubes as necessary

Adjust Volume and Speed appropriate to cell concentration².

Adjust stopping gate and Record samples

Cleaning

Run Sanitize SIP from *Instrument Tab*. Choose the appropriate option if using plates

Wait until cleaning has completed and rinse.

LAST USER: Run Shutdown from *Instrument Tab*³.

Exit software then log off computer. End Kiosk session.

¹If another user is logged in, restart computer. Ensure no plates are running.

²Evs/sec<35K. For accurate cell counts, evts/sec<8K.

³When progress bar reaches ~6%, you can log off the computer. DO NOT SHUTDOWN computer.