
USING THE PLATE HOLDER ON THE FACSMELODY


UNC FLOW CYTOMETRY CORE FACILITY

EJECT PLATE HOLDER


1. Insert **ACDU** into Sort block



2. Navigate to “2 Set Up Sort” > Select Plate and define sorting parameters

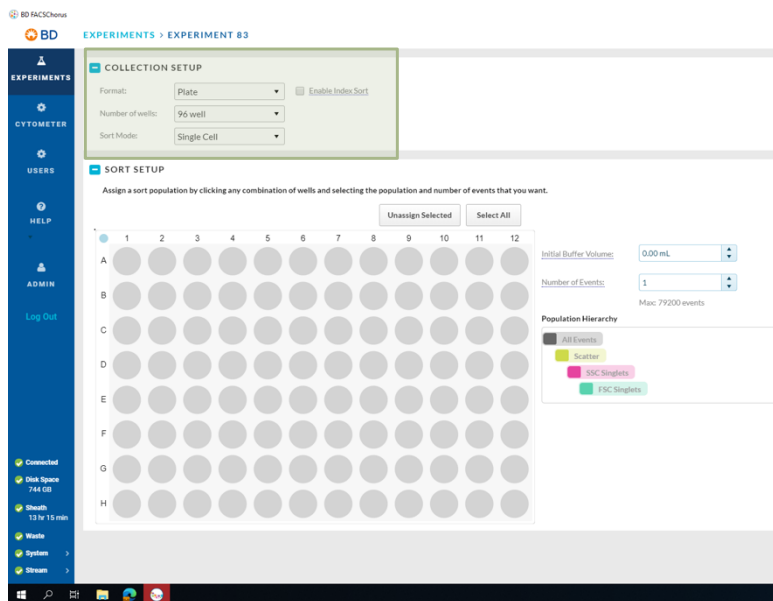

3. Navigate to “3 Sort” >Select **EJECT** in the Sort Status Menu

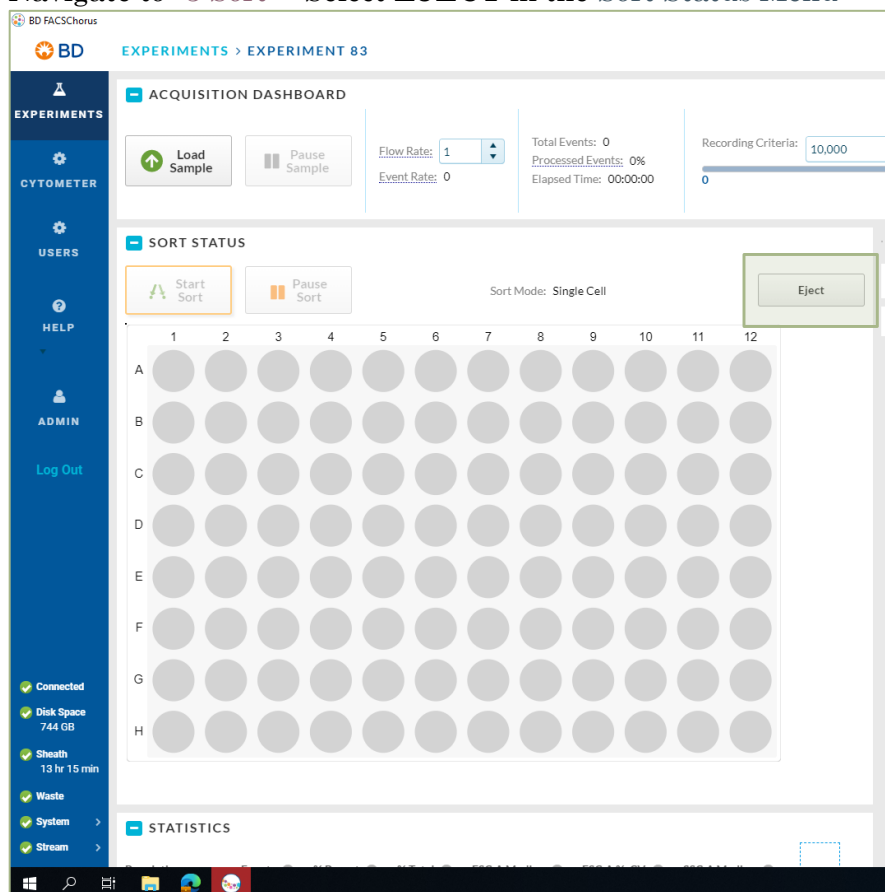


4. Place Plate on Plate holder and proceed to sort

RETRACT PLATE HOLDER

1. Keep **ACDU** in Sort Block



2. Navigate to the “2 Set up Sort” > Select TUBES

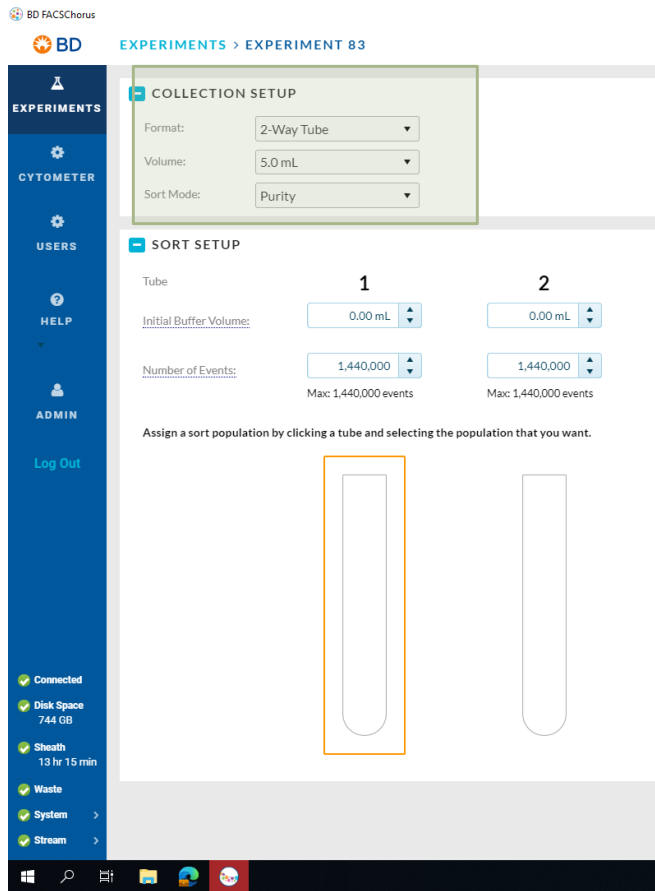

3. Navigate to “3 SORT” > Select **RETRACT** in Sort Status Menu

